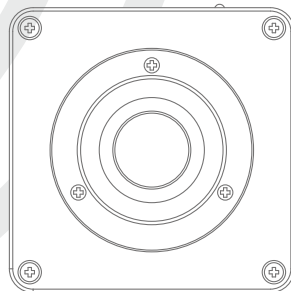
Moticam_i QUICK START GUIDE

EN Quick Start Guide

Thank you for purchasing this Moticam. This camera is a delicate instrument and should be handled with care. When not in use, please store the camera in a safe place with the dust cap attached. Do not let the camera get in contact with water or expose it to extreme temperatures (below 0 or above 50 degrees Celsius). In order to receive the highest level of service, please retain your receipt as proof of purchase and **register your camera online at www.motic.com today.** Should you have any questions please do not hesitate to **contact our customer service.**

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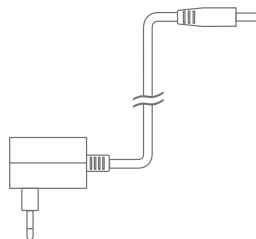
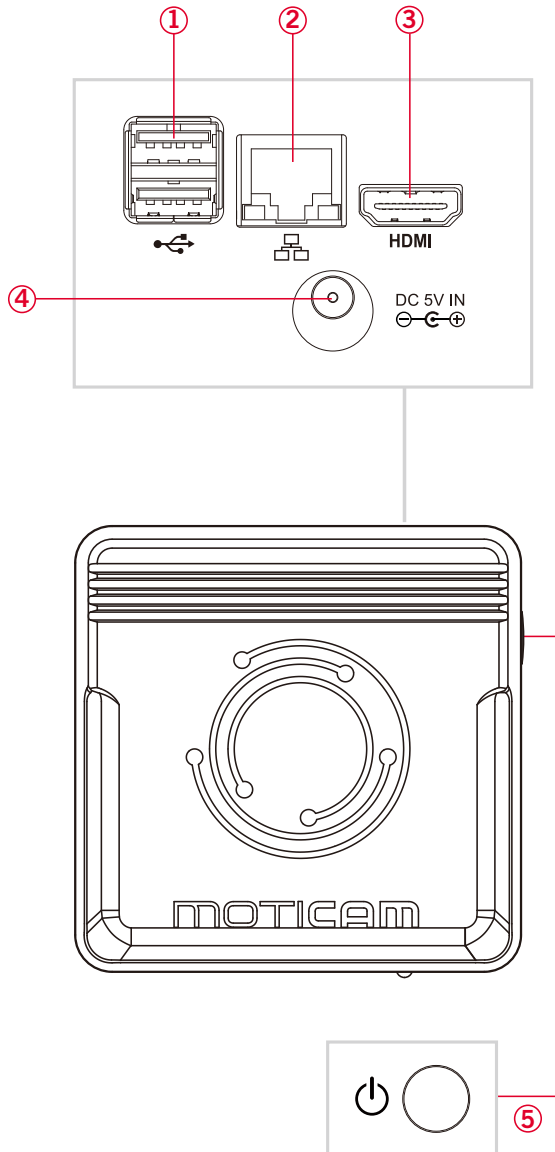
EN Controls and Ports on the Camera

1. The USB connection can be used to attach a remote keyboard and mouse or a memory stick to the camera. It can also be used to connect an optional Wi-Fi antenna.
2. The RJ-45 port can be used to network the camera with a router, it can also be used to connect the camera to a computer running Motic Images Plus.
3. The HDMI port can be used to connect to an HDMI monitor. A cable is included in the package.
4. Please use only the supplied power adapter to power the Moticam I.
5. The on/off button controls the power to the camera. When the camera is on, an indicator light will illuminate on the front of the camera.

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EN On-Screen Display

When the camera is connected to an HDMI display and controlled through a USB mouse, you can simply move the mouse cursor to the right to activate the on-screen function display for capturing images and also for adjusting various aspects of the image parameters.

You can also use this setting to change the on-screen display language from the default English to Japanese or Chinese.

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EN Using the camera on a Trinocular Microscope

1. Install the Motic Images Plus application program from the CD included.
2. Remove the Dust Cap from the camera (A) and attach the microscope's C-Mount adapter (*Please note that the C-Mount adapter is specific to your microscope. For Motic microscopes, please consult the relevant Motic Accessory list and choose an appropriate adapter for your Moticam. Depending on the type of adapter and microscope, the C-Ring (B) may be necessary between the camera and the C-mount adapter).
3. Connect the HDMI cable (E) into an HDMI screen and you can also connect an optional remote mouse and keyboard into one of the USB ports. The other USB port will accept a USB memory stick. You can connect the camera to a computer through the RJ-45 port and run it with the supplied Motic Images software.

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